



Title: *Hybridoma Media Preparation: DMEM-10*

No: RTLP-GL-Ab-6

Location:
Old CCRC Tripp Lab

Approval Date:
10 September 2004

Supersedes Date:

Materials:

Page 1 of 1

- | | | | |
|----------------------------------|----------------------------|----------------------|------------------|
| •Lab coat | •Fetal Bovine Serum (FBS) | •DMEM | •Pipettes |
| •Gloves | •Sterile laminar flow hood | • Pen/Strep Solution | •Pipetteman |
| •Gentamycin | | •L-Glutamine 200mM | •Pipette Aid |
| •Phosphate Buffered Saline (PBS) | | | •Pipetteman tips |

Procedure:

DMEM-10:

- Under sterile conditions combine 1.0 mL Gentamycin, 10 mL Pen/Strep, and 100 mL Hyclone FBS to 1L DMEM.
- Date and initial the tube and label the contents.
- Store at 4°C

NOTE: L-Glutamine is no longer effective after a two week time period, so use accordingly.

Author	Management Approval/Date	Effective Date