



Title: *Hybridoma Media Preparation: 50X HAT
Supplement*

No: RTLP-GL-Ab-8

Location:
Old CCRC Tripp Lab

Approval Date:
10 September 2004

Supersedes Date:

Materials:

Page 1 of 1

- | | | | |
|-----------------|------------------|------------------|----------------|
| •Lab coat | • 10 mL syringe | •Sterile laminar | •Pipettes |
| •Gloves | •18 gauge needle | flow hood | •Pipette Aid |
| •HAT supplement | •5 mL tubes | | •-20°C Storage |
| •DMEM | | | |

Procedure:

- Using a 10 mL syringe and 18 gauge needle, add 10 mL of sterile DMEM to one bottle of HAT supplement.
- Invert bottle several times to mix into solution.
- Remove the metal criped rim and aseptically remove the top of the HAT supplement.
- Aliquot in 2 mL amounts into sterile tubes.
- Date the tube and label the contents.
- Store at -20°C.

Author	Management Approval/Date	Effective Date

