



**Title:** *Buffers and Stock Solutions: NaOH, 10M*

No: RTLP-GL-BSS-9

Location:  
*Old CCRC Tripp Lab*

Approval Date:  
10 September 2004

Supersedes Date:

**Materials:**

Page 1 of 1

- Lab coat
- Gloves

- NaOH
- dH<sub>2</sub>O

- Analytical balance
- 100 mL beaker

- Stir plate
- Stir bar

**Procedure:**

1. Weigh out 400 grams NaOH and bring up to solution with 1L dH<sub>2</sub>O.
2. Stir to mix into solution.

Author	Management Approval/Date	Effective Date